
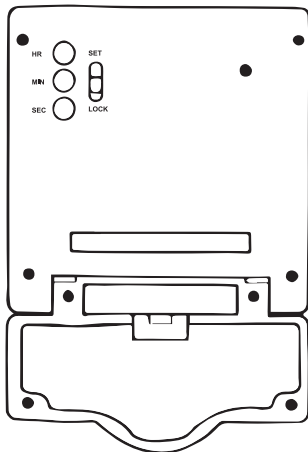


# **White Large Display Timer**

**810005  
Instruction Manual**

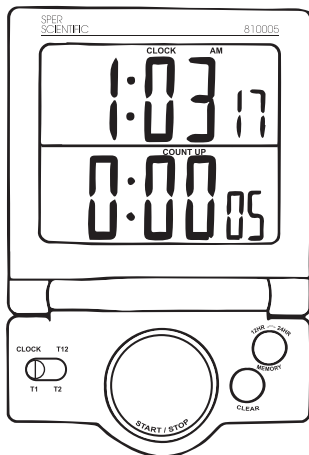

## OPERATING INSTRUCTIONS

### To Set Clock:

12HR (a.m./p.m.) 24HR International time display format user selectable. Slide switch to **CLOCK**, press **12HR** ⇌ **24HR** button for desired format. Depress **HR**, **MIN**, **SEC** buttons at the unit back to actual time when the slide switch is in **SET** mode. After setting, slide to **LOCK** mode to avoid unintentional touch of HR, MIN, SEC buttons resulting in wrong readouts.

## Count-Up:

When the front sliding switch is in CLOCK mode, press START/STOP once to start, second press to pause, third press to restart. Press CLEAR after pausing will reset digit to 00:00:00. Count-up to 20 HR... Maximum count-up time capacity is 19:59:59.

## Timer 1 & 2

Slide front switch to **Timer 1** or **Timer 2**, upper or lower row will show 0:00:00. When flashing depress **HR, MIN, SEC** buttons at the back of the unit to actual timer when slide switch is in **SET** mode, after Setting, slide to LOCK mode to avoid unintentional touch of HR. MIN, SEC buttons.

## Memory

Slide front switch to **Timer 1** or **Timer 2**, upper or lower row will show 0:00:00. When flashing depress **HR, MIN, SEC** buttons at the back of the unit to actual timer when slide switch is in **SET** mode, then press **MEMORY** switch, MEMORY icon is on. After setting, slide to **LOCK** mode to avoid unintentional touch of HR, MIN, SEC buttons.

## **WARRANTY**

Sper Scientific warrants this product against defects in materials and workmanship for a period of **five (5) years** from the date of purchase, and agrees to repair or replace any defective unit without charge. If your model has since been discontinued, an equivalent Sper Scientific product will be substituted if available. This warranty does not cover, batteries, battery leakage, or damage resulting from accident, tampering, misuse, or abuse of the product. Opening the meter to expose its electronics will break the waterproof seal and void the warranty.

To obtain warranty service, ship the unit postage prepaid to:

### **SPER SCIENTIFIC LTD.**

8281 East Evans Road, Suite #103  
Scottsdale, AZ 85260

The defective unit must be accompanied by a description of the problem and your return address. Register your product online at [www.sperwarranty.com](http://www.sperwarranty.com) within 10 days of purchase.